**Bürkle GmbH**

Rheinauen 5 | D - 79415 Bad Bellingen  
Telefon +49 (0) 76 35 8 27 95-0  
Telefax +49 (0) 76 35 8 27 95-31  
info@buerkle.de | www.buerkle.de

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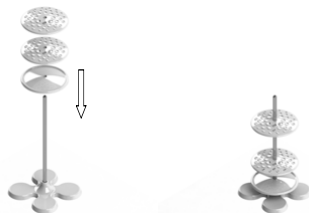
## General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

## Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Use only in ventilated rooms or under a suitable exhaust hood.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ The instrument is suitable for indoor use.
- ▶ The instrument is designed for stationary use and is not suitable for the transport of pipettes.

## Safety instructions during use

### **WARNING!**

Danger due to unsuitable working area. Danger of tipping over.

- ▶ Bright, dry and clean working place at table height without obstructions.
- ▶ Horizontal stand of the frame with sufficient space.
- ▶ The instrument has to be mounted on a horizontal, even surface.
- ▶ The device can be used in the temperature range of 0 - 40 °C.
- ▶ The instrument is suitable for the use with pipettes, thermometers and test tubes.

## Layout and function

- ▶ The instrument consists of a foot, a drip pan, two perforated plates made of polypropylen (PP) and a support rod made of chrome-plated steel.
- ▶ The instrument is suitable to hold pipettes, thermometers and test tubes. It serves for storing, draining and drying pipettes.

- ▶ The instrument is suitable for 40 pipettes with a maximum diameter of 18 mm. Number and diameter of the openings:
  - 4 x Ø 18 mm
  - 4 x Ø 16 mm
  - 8 x Ø 14 mm
  - 8 x Ø 12 mm
  - 16 x Ø 10 mm
- ▶ The instrument is maintenance-free.

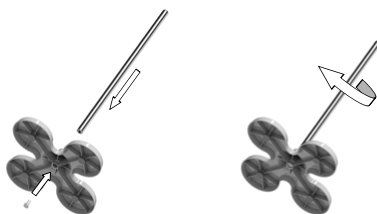
## Scope of delivery

1 x	Support rod, Ø 16 mm, chrome-plated steel	①
1 x	Foot, PP	②
1 x	Drip pan with magnets, PP	③
2 x	Perforated plates with magnets, PP	④
1 x	Directions for use	

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

## Assembly

1. Screw off the screw from the support rod ①.
2. Insert the support rod from above into the foot ②.
3. Fasten the support rod with the aid of the screw from below (by rotating the support rod).

4. Plug the drip pan ③ onto the support rod and slide it downwards.
5. Plug the perforated plates ④ onto the support rod and move them to the desired height.

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